100% Fiocchetti the new standard

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- + 50 years in the refrigeration field to reach "the cold management at the state of the art"
- + Main sectors: medical, pharmacies, hospitals, blood transfusion centers
- + Innovative, functional design
- + Latest generation electronics, highest control on the performance of the equipment and on power consumption
- + Highly professional refrigerators, worldwide present thanks to our distributors

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The professional refrigerators range MEDIKA 2T have been designed specifically for pharmacies and pharmaceutical sector: two temperature settings, two cooling units and two totally independent chambers to store at best different groups of products. Each chamber can be set independently between +2° and +15°C and is provided as a standard with acoustic and visual alarms. This range offers units with different capacities, from 280 to 1500 liters, all supplied with triple layer glass doors and can be customized as far as the internal fittings (with shelves and/or telescopic drawers). The range MEDIKA 2T complies with Italian pharmacopeia and European drugs storage normative and it's available in ECT-F, ECT-F PLUS or ECT-F TOUCH version.

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B+B= Lamiera plastificata bianca antibatterica interno ed esterno / Pre-coated antibacterial sheet steel, white colour, both inside and outside

B+X= Lamiera plastificata bianca antibatterica esterno e acciaio inox AISI 304 interno / Pre-coated antibacterial sheet steel, white colour, outside and AISI 304 stainless steel inside

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Sonda umidità
Humidity probe

To complete an already rich standard equipment, the series MEDIKA, MEDIKA 2T and LABOR can be customized with the following accessories:

- + **Wire shelf** in a-toxic plastic coated steel, supported by anti-tilt clips.
- + **Aluminium drawer**, totally extractible on telescopic slides, adjustable in height, bottom made of painted steel, bee-hive structure type, supplied with adjustable dividers made of polypropylene. The front is fitted with an ergonomic aluminium handle.
- + **Stainless steel wire shelf**
- + **Stainless steel wire drawer** on telescopic guides
- + **Drawer** for narcotics with keylock
- + **Stainless steel basket** mounted on telescopic slides.
- + **Stainless steel rack**, dimensions cm. 14x42/50x15
- + **Castors kit**, nr.2 with brakes
- + **Electric digital key lock**: the door opening is possible by pressing the icon "key" on the touch screen. This function can be password protected to avoid access from un-authorized personnel. In case of power failure the door can be opened manually using a supplied tool
- + **External terminal board** for dry contact, Ethernet and RS485 ports
- + **Back up batteries kit** for acoustic and visual power failure alarm: this accessory includes nr 8 eco-friendly batteries Ni-MN type, easily replaceable from end users
- + **DMLP Digital Monitor Touch**: independent plug-in module for temperature recordings and alarms on SD card even during power failure, working with PT 100 class A probes and with a 30 seconds sampling. This module includes nr 2 dry contacts to remote alarms and a back-up batteries kit. The DMLP compares continuously the data coming from its PT100 probes and those of the ECT-F Touch controller, signaling any eventual discrepancy. This check allows to be informed immediately if any of the thermoregulation probes are unbalanced and thus to grant the perfect storage of products.

- + **Wi-Fi Module**: plug-in accessory for monitoring, connection and control at the distance of the equipment
- + **Gsm Communicator**: plug-in module for the notification, via SMS, of the functioning and possible faulty condition of the equipment. Every ECT-F Touch controller can be equipped with a GSM module, thus becoming an independent unit sending and receiving different sort of SMS. Monitoring SMS: calling the SIM card of the refrigerator will generate an SMS containing all functional data. ALERT SMS: failures of faulty conditions are notified to the pre-set numbers. When equipped with back-up batteries, in case of power failure, the GSM will inform of this condition and of power restore, once this will be the case. COMMAND SMS: messages can be sent to modify the controller parameters, the turning on or off of the unit, to change the set-point
- + **Voltage stabilizer**
- + **DIN 58345 conformity kit**: along with the equipment standard features (for instance: safety thermostat), it included the back-up batteries kit, electric digital key lock, the external terminal board for dry contact, Ethernet and RS485 ports.
- + **Humidity probe**: to monitor and record the humidity level values inside the compartment
- + **Temperature product simulation** (ballasted probe) - PT 100 probe
- + **PT100 probe** (free contacts)
- + **Pass-through port with cap**, diam. mm 35 (as a standard, placed on right side)
- + **Pass-through port** mm 15 (as a standard, placed on the back side)
- + **Internal socket** 220V
- + **Double internal socket** 220V with ON/OFF switch
- + **Temperatures certificate**: test report, minimum nr.9 probes for 12 hours
- + **Wooden crate package**

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